

CI-muINTEPI

Cat. No.: INS-CI-1007



Biosafety Level
Level 1



Storage
vapor phase of liquid nitrogen



Culture Conditions
muINTEPI Medium (INS-ME-1005)
Collagen solution (INS-SU-1017)

General Information

CI-muINTEPI is a murine intestinal epithelial cell line from mouse E19.5 fetuses. The cell line was immortalized using the [CI-SCREEN](#) technology ([Lipps et al. 2018; Nat Comm](#)) and characterized in [Schwerk et al. 2013](#).

Organism: Mus musculus (C57BL/6, "Black6")

Tissue: intestine

Growth properties: adherent

Cell culture media and reagents

Product	Cat. No.	Volume
muINTEPI Medium (includes basal medium and supplements)	INS-ME-1005	500ml
Collagen solution	INS-SU-1017-50ml INS-SU-1017-300ml	50ml 300ml
Freezing medium	INS-SU-1004	30ml

Note: The medium does not contain antibiotics. However, it may simply be supplemented with standard antibiotics.

Intended Use

This product is intended for in vitro research use only. It is not intended for any animal or human therapeutic or diagnostic use.

Quality control

Each vial contains $\geq 5 \times 10^5$ cells. Viability is $\geq 80\%$. Cells are negative for mycoplasma contamination.

Upon arrival

Cells are routinely shipped on dry ice. Check all containers for leakage and breakage. Check if cells arrived frozen. After arrival, store the cryopreserved cells in liquid nitrogen vapor, or seed them immediately (please see page 2).

Note: Cells may be stored at -80°C for short periods (<2 days), but this results in reduced viability and irreversible cell damage.

Medium storage and preparation

Medium including supplements is shipped cooled at $4-8^\circ\text{C}$. Store reagents according to the instructions below upon arrival.

Material:

- muINTEPI Medium (INS-ME-1005). Includes Basal Medium (500ml) and Supplements (30ml; INS-ME-1005BS).

Storage:

- Store Basal Medium at $4-8^\circ\text{C}$
- Store Supplements at -20°C
- Store completed medium (Basal Medium plus Supplements) at $4-8^\circ\text{C}$. Completed medium is stable for at least 1 month at $4-8^\circ\text{C}$.

Protocol:

- 1) Thaw Supplements at $15-25^\circ\text{C}$.
- 2) Add 30ml Supplements to 500ml Basal Medium and store at $4-8^\circ\text{C}$. Completed medium is stable for at least 1 month at $4-8^\circ\text{C}$.

Note: Supplements may be aliquoted and stored at -20°C before completing the medium. For example, aliquot $5 \times 6\text{ml}$ and then add 6ml Supplements to 100ml Basal Medium.

Collagen coating of cell culture vessels

Material:

- Cell culture vessel(s)
- PBS
- Collagen solution (INS-SU-1017)

Protocol:

- 1) Cover the cell culture vessel with coating solution (see table below for the required volume).
- 2) Incubate the cell culture dish for at least 2h (up to overnight) at 37°C in the incubator.
- 3) Aspirate coating solution.
- 4) Wash once with PBS and aspirate PBS.

Vessel	Surface area (cm^2)	Volume (ml)
T75	75	2.5
T25	25	1.4
6well	9.6	0.7
12well	3.5	0.25
24 well	1.9	0.1
96well	0.32	0.05

Note: Coated cell culture vessels may be stored sealed at $2-8^{\circ}\text{C}$ for up to 7 days.

Recover cryopreserved cells

Do not thaw the cells until the recommended medium and coated flasks are on hand. For initial recovery (after delivery), we recommend thawing the cells on a T25 flask.

Material:

- complete medium
- 15ml tube
- Collagen-coated culture vessel

Protocol:

- 1) Add 4ml pre-warmed medium to a 15ml tube.
- 2) Quickly thaw the cryovial at 37°C in a water bath until only a few ice crystals are visible. Disinfect vial briefly with 70% Ethanol.

- 3) Transfer thawed cell suspension to the 15ml tube containing 4ml medium. Avoid excessive pipetting up and down.
- 4) Centrifuge cells at $200 \times g$ for 4min.
- 5) Aspirate supernatant.
- 6) Gently resuspend the cell pellet in complete medium.
- 7) Transfer cells in coated cell culture vessel and place in the incubator (37°C , 5% CO_2).
- 8) Change the medium after 2 days.

Routine Subculture

Change medium every 2-3 days and split the cells at 70-90% confluence. The split ratio after recovery from cryopreservation should not exceed 1:2. For routine maintenance, split ratio can be increased to 1:5 to 1:10.

Material:

- complete medium
- PBS
- Trypsin/EDTA solution (TE)
- Collagen-coated culture vessel

Protocol:

- 1) Aspirate medium.
- 2) Wash with PBS and aspirate PBS.
- 3) Add Trypsin/EDTA (TE) solution to the cells and incubate at room temperature or 37°C for 5-10min, or until the cells detach.
- 4) Examine the cells under a microscope. When the cells start to detach, gently tap the side of the vessel to loosen the remaining cells.
- 5) Resuspend cells in complete medium thereby inactivating the Trypsin/EDTA (TE) solution.
- 6) Transfer an aliquot of the cell suspension to a new coated cell culture vessel containing fresh complete medium.
- 7) Incubate at 37°C and 5% CO_2 .

Vessel	Medium or PBS (ml)	TE (ml)
T75 flask	8-10	2
T25 flask	4-5	1
6well	1.5-2	0.5
12well	1	0.2
24 well	0.5	0.1
96well	0.1	0.05

Cryopreservation

Cell should be grown to 90% confluence before cryopreservation. Avoid full confluence before cryopreservation.

Material:

- Freezing medium (INS-SU-1004)
- PBS
- Trypsin/EDTA solution (TE)
- 2% FBS in PBS
- 15ml tube
- cryovial(s)
- freezing container ("Mr. Frosty" or similar)

Protocol:

- 1) Aspirate medium.
- 2) Wash with PBS and aspirate PBS.
- 3) Add Trypsin/EDTA (TE) solution to the cells and incubate at room temperature or 37°C for 5-10min, or until the cells attach.
- 4) Examine the cells under a microscope. When the cells start to detach, gently tap the side of the vessel to loosen the remaining cells.
- 5) Resuspend cells in 2% FBS in PBS and transfer to a 15ml tube.
- 6) Centrifuge cells at 200×g for 4min.
- 7) Aspirate supernatant and gently resuspend cell pellet in Freezing medium (approx. 1 × 10⁶ cells/ml).
- 8) Transfer cell suspension into cryovial(s) and place them into a freezing container ("Mr. Frosty" or similar).
- 9) Place the freezing container at -70 to -80°C for 16-24h.
- 10) Transfer cryovials to liquid nitrogen vapor for long-term-storage.

Contact Information

InSCREENeX GmbH
 Inhoffenstr. 7
 38124 Braunschweig
 Germany

Email: info@inscreenex.com

Website: inscreenex.com

Phone: +49 531 6181 5080

Fax: +49 531 6181 5002

Selected References

Cell line generation and characterization:

Schwerk J et al. Generation of mouse small intestinal epithelial cell lines that allow the analysis of specific innate immune functions. *PLoS One*. 2013;8(8):e72700. Published 2013 Aug 5. doi:10.1371/journal.pone.0072700

Heterogenous type-III interferon response:

Bhushal S et al. Cell Polarization and Epigenetic Status Shape the Heterogeneous Response to Type III Interferons in Intestinal Epithelial Cells. *Front Immunol*. 2017;8:671. Published 2017 Jun 12. doi:10.3389/fimmu.2017.00671

Interferon-λ-induced transcriptional profile:

Selvakumar TA et al. Identification of a Predominantly Interferon-λ-Induced Transcriptional Profile in Murine Intestinal Epithelial Cells. *Front Immunol*. 2017;8:1302. Published 2017 Oct 16. doi:10.3389/fimmu.2017.01302

Cryptosporidium Infection:

Li M et al. Induction of a Long Noncoding RNA Transcript, NR_045064, Promotes Defense Gene Transcription and Facilitates Intestinal Epithelial Cell Responses against Cryptosporidium Infection. *J Immunol*. 2018;201(12):3630-3640. doi:10.4049/jimmunol.1800566

Complement system:

Sünderhauf A et al. Regulation of epithelial cell expressed C3 in the intestine - Relevance for the pathophysiology of inflammatory bowel disease?. *Mol Immunol*. 2017;90:227-238. doi:10.1016/j.molimm.2017.08.003